Shodex

Operation Manual

(temporary)

Anion separation columns for suppressor method Shodex IC SI-50 4E

Showa Denko K.K.

Introduction

The Shodex IC SI-50 4E column is developed for use in suppressor method anion chromatography.

The seven anions (fluoride, chloride, nitrite, bromide, nitrate, phosphate and sulfate) and organic acids (for examples: acetate, formatecan, methacrylate and oxalate) be separated with high sensitivity.

Specifications

Table 1

Product name	Column size	Theoretical plate	Particle	Application
	(i.d. x L (mm))	number	size	
IC SI-50 4E	4.0 x 250	≥10,000 (Br)	5 um	For analysis
IC SI-90G	4.6 x 10	****	9 um	As guard column

The packing material is an anion exchanger made from polyvinylalcohol gel chemically bonded with quaternary ammonium.

The liquid with which the columns are filled at delivery is a mixture of 3.2 mM Na_2CO_3 and 1.0 mM $NaHCO_3$.

Table 2

	SI-50 4E
Column material	PEEK
Recommended eluent	3.2 mM Na ₂ CO ₃ + 1.0 mM NaHCO ₃
Maximum flow rate	0.8 mL/min
Recommended flow rate	0.7 mL/min
Maximum pressure	15.0 MPa
Usable pH range	pH 3~12
Recommended temperature range	20 ~ 60 °C

Sample pretreatment

- 1) Inject the sample into the column only after it has been passed through a 0.45 μ m membrane filter to remove particles.
- 2) Any sample containing protein should be injected into the column only after protein has been eliminated from the sample.
- 3) Inject the sample containing organic impurities into the column only after the sample has undergone solid extraction treatment (Sep-Pak PS-1).

Eluent

Normally aqueous solution of descrived in the tabel 2 can be used as an eluent for Shodex IC SI columns.

1) 3.2 mM $\rm Na_2CO_3$ + 1.0 mM $\rm NaHCO_3$ Measure 0.339g of $\rm Na_2CO_3$ and 0.0840g of $\rm NaHCO_3$ into a 1 liter measuring flask. Make it up a 1 liter solution using distilled and deionized water.

Storage

The column should be throughly flushed with fresh eluent. Column disconnected from the LC system should be tightly capped both ends to prevent internal drying, and stored in an room that has less temperature fluctuation.

Regenaration

Cause	Washing procedure	
	Washed by the following steps. (flow rate 0.3 ml/min)	
Polution by low valency	1. 25 minutes : deionized water	
hydrophilic ions	2. 100 minutes: 10 times concentrated eluent	
	3. 25 minute: deionized water	
	4. 100 minutes: eluent	
Washed by the following steps. (flow rate 0.3 ml/mi		
Polution by high	1. 25 minutes : deionized water	
valency hydrophobic	2. 20 minutes: 5% acetonitrile	
ions	3. 100 minute: 100% acetonitrile	
	4. 50 minutes: deionized water	
	5. 100 minutes: eluent	