# Operation Manual

## Shodex IC YS-50

L-25

(Please read this manual carefully before using the column to keep its performance and life on a standard level.)

### 1. Introduction

Shodex IC YS-50 is an ion chromatography column suitable for simultaneous analysis of univalent and divalent cations. It has the merit of being able to detect a trace ammonium ion under a high concentration of sodium ion. It is used for both suppressor method and non-suppressor method.

### 2. Instructions in handling <Important>

#### Caution!

\* Take notice of keeping instructions about the solvents and the reagents used with the column not to occur problems related to losing your health or leaking.

#### Attention!

\* Use the column within the regular range of flow rate, pressure and temperature. There is a danger of deteriorating the performance when it is handled beyond the permissible range even for a short time. See the clause "Usable conditions" about the permissible range.

### 3. Specifications

Column size	: 4.6 mm I.D. × 125 mm L.
Column material (inside)	: SUS 316
Packing material	: Porous particle of poly(vinyl alcohol) type bonded with carboxyl group
In-column solvent (initial)	: 4.0mM methane sulfonic acid
Number of theoretical plates	: $\geq 5,500$ per column ( $\text{Ca}^{2+}$ )

### 4. Usable conditions

Flow rate	: $\leq 2.0$ ml/min
	* It should be 1.0ml/min or less for replacing the in-column solvent.
Pressure	: $\leq 10$ MPa per column
Temperature	: 20~60 °C
pH	: 2~12
Eluent	: 4.0mM methane sulfonic acid is usually used.
	* Acetonitrile can be added within 30 vol%.

#### Attention!

- 1) Replace the solvent in the chromatograph with the eluent to be used before connecting the column.
- 2) Connect the column so that the flow direction corresponds to the arrow mark on the tag.
- 3) Connect the column with a male screw and a ferrule made of resin. Pay attention not to screw too tightly. PEEK connector (a ferrule unified with a male screw) which needs no tool is recommended.
- 4) Filtrate the sample with a disposable filter (0.45  $\mu\text{m}$ ) to prevent deterioration by adsorbing insoluble matters.
- 5) Install a guard column (YS-G) upstream to protect the main column from deterioration with adsorbates.
- 6) When the column is not used for a month or more, replace the in-column solvent with the standard eluent, close each end with a stopper, and store it at room temperature.
- 7) Do not remove the end fittings of the column under any circumstances.
- 8) Do not make a strong impact on the column: such as hitting or dropping on the floor.

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